

Heat Inactivation of FBS

Materials:

500 mL bottle of FBS (VWR/Seradigm or other current vendor)

1. Incubate in a water bath at room temperature until thawed (~2h).
2. Turn on water bath with a set point of 56°C.
3. Once set point is reached, incubate in 56°C water bath:
 - a. 50 mL for 25 min
 - b. 500 mL for 45 min
4. Transfer to 4°C for ~30 min to chill.
5. Aliquot into 50 mL conical tubes labeled with: initials, product #, lot #, date. Use laser printed cyro babies tags if desired.